


Head-mounted condenser microphone and power adapter

- There are two leads from the microphone. The light grey plug with the microphone symbol should be connected to the microphone power adapter input (not the black plug). The microphone power adapter needs to be switched on (thumb wheel on the side of the adapter).
- The microphone power adapter output should then be switched on and connected to the microphone input on the EPG3 Scanner unit.
- The microphone gain knob on the front of the SPI unit should be around the 10 o'clock mark.
- Then you should see movement on the audio VU meter on the front of the SPI unit.

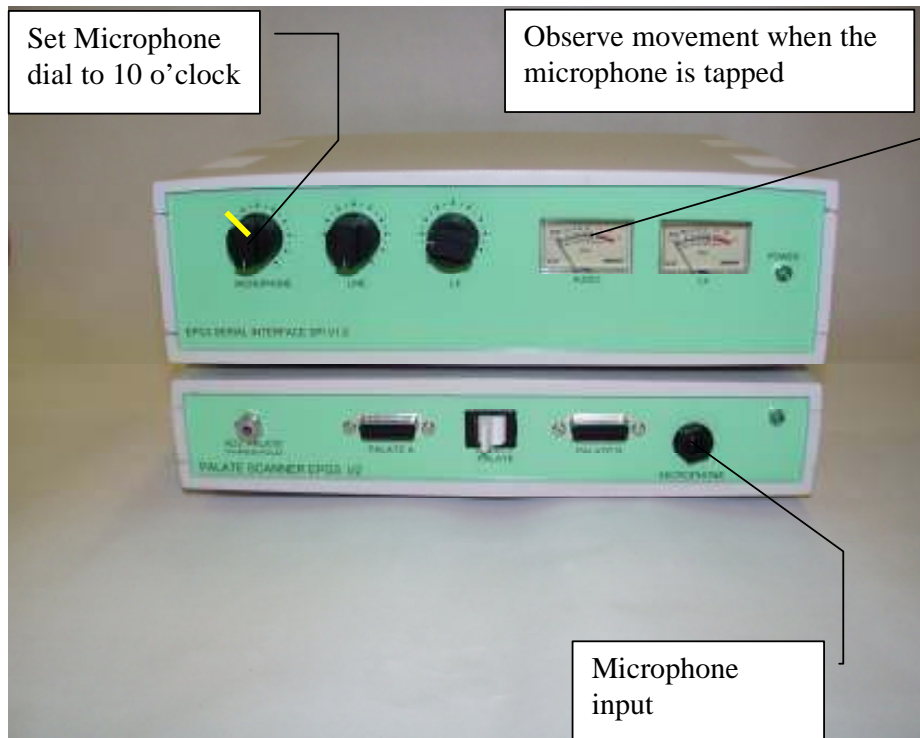

- Connect the “Audio out” and “Lx out” phono sockets to the “Line-in” socket (usually blue) on the PC soundcard with the twin phono to minijack cable provided.



- Go to the Options|Audio menu to select the audio mixer and set the Recording

